

MESOTRON

Quartz table stopclock

zu beziehen bei

sold by

www.conatex.com



Preparation

Remove the battery-cover located under the base. When changing the battery type AA of 1,5 V, respect polarity. Replace battery cover. Reset the stopwatch by pressing the black button.

Functions

1. start: press the red button
stop: press the red button
reset: press the black button
2. addition: press the red button consecutively as often as required
3. tempo: start-reset with automatic starting of the second hand. Press red button and then the black one as often as requested.

Technical data

| | |
|-------------------------|---|
| movement: | quartz controlled 32,768 Khz |
| accuracy: | +/- 14 sec./month |
| dial: | first hand 60 min. second hand 60 sec. - 100/100 min. diameter 110 mm ø |
| time units: | 1 sec. and 1/100 min. |
| case: | ABS, 175 x 130 x 40/95 mm |
| weight: | 370 g approx. |
| battery: | Mignoncell 1,5 V (AA) |
| duration: | 1 1/2 years |
| battery life indicator: | in case of critical battery capacity, the hand is not running smoothly (0,75-1,25-0,75-1,25...) sec. Then there is only power left for about 1 week |
| temperature in use: | -10° C to + 60° C |

This product is German made.